**Compact Dry CF medium for coliforms/ Medium für Coliforme / milieu pour coliformes / medio para coliformes / medio per coliformi / meio para coliformes**

40 plates/Platten/plaques/placas/ lastre/placas	ID-No. 1 000 867
240 plates/Platten/plaques/placas/lastre /placas	ID-No. 1 000 868
500 plates/Platten/plaques/placas/lastre /placas	ID-No. 1 400 868
1400 plates/Platten/plaques/placas/lastre /placas	ID-No. 1 502 881

**English****Deutsch****Français**

*Compact Dry CF is a ready to use, chromogenic plate for the enumeration of coliforms*

**Sample pretreatment****Viable count in water or liquid foodstuff**

Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

**Viable count in solid foodstuff**

Add buffer solution (e.g. 0.9 % saline solution, PBS, buffered peptone water, etc.) to the sample and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) on the middle of the dry sheet of the Compact Dry plate.

**Viable count in swab test specimen**

Use the swab to wipe the surface, put into the device with wiping solution. After shaking the device (approx. 10 sec) drop 1 ml of wiping solution (dilute if necessary) on the middle of the Compact Dry plate. It is recommended to use "Swab for Compact Dry" offered by HyServe Id-No. 1 002 952/3 (40/240 pieces)

**Test instructions**

1. Open the cap and drop 1 ml of specimen on the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. Put the cap again on the plate and write the information needed on the memorandum section.
4. Turn over the capped plate and put in the incubator.
5. After incubation count the number of colored colonies underneath the plate. White paper placed under the plate helps to count the colonies.

**Incubation time** 24 ± 2 hours

**Incubation temperature**

**Microval** 37 ± 1°C

**AOAC** 35 ± 2°C

**Interpretation of the results**

Colonies grown are blue/blue green colored.

**Storage and shelf life**

Keep at room temperature (+ 1 to + 30 °C). Total shelf life 24 months after manufacturing.

**Notes**

- The growth of bacteria other than coliforms is mainly inhibited.
- After use please follow the current disposal regulations.
- High concentrations on plates (> 250 CFU) will cause the entire growth area to become white/blueish. In this case dilute the sample.
- The growth area is 20 cm<sup>2</sup>. The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, the bacteria count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids. However, dilutions are recommended.
- Compact Dry plates are produced at an ISO 9001/ISO 13485: 2003 certified site.
- AOAC approved; certificate No. 110401
- MicroVal approval No.0806-003LR/ISO 4832 (2006)
- ISO EN 16140:2003
- NordVal approval No. 035

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